



ThermoFisher
S C I E N T I F I C

Contamination Analysis and Identification by Vibrational Spectroscopy

Matt Bartucci

2017 NASA Contamination, Coatings, Materials, and Planetary Protection Workshop

NASA Goddard Space Flight Center

The world leader in serving science


Deformulation and Contamination Analysis via FTIR

- Reverse Engineering
- Failure Analysis
- Material Analysis

- Polymers
 - Plastics
 - Fillers
 - Rubbers (!)
 - Carbon Black: O-Rings

- Epoxies, Resins, Adhesives

- Excellent 1st approach technique

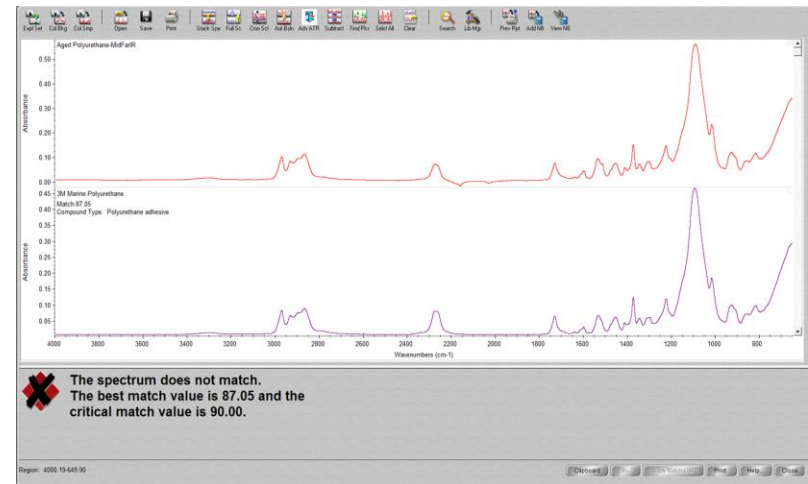

- **ATR-IR**

- Contamination analysis of organic and inorganic materials
- Reverse engineering of heat shrink tubing



- **TGA-IR**



- Failure analysis of o-rings




Basics: Depth of Penetration

- Attenuation occurs by sample absorption: $A = \epsilon lc$
- The Evanescent wave probes short distance (~1-2 micron)
 - Measures surface & near-surface only
 - Path Length 'l' characterized by depth of penetration

$$d_p = \frac{\lambda}{2\pi n_{\text{atr}} \left[(\sin^2 \theta) - \left(\frac{n_{\text{sample}}}{n_{\text{atr}}} \right)^2 \right]^{1/2}}$$


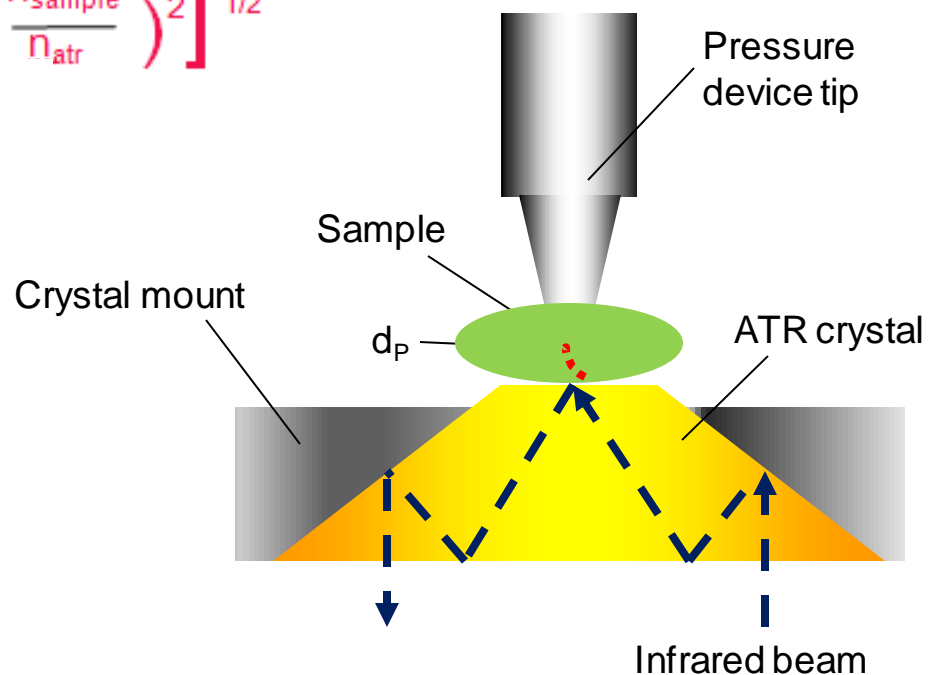
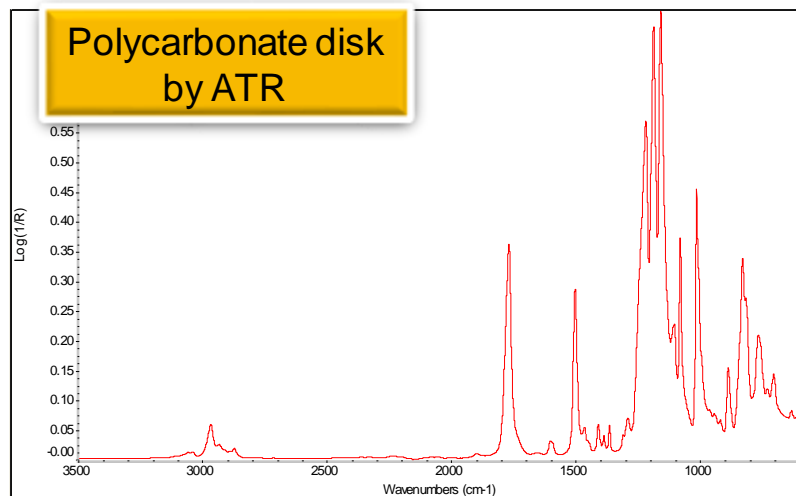


Depth of Penetration with Wavelength

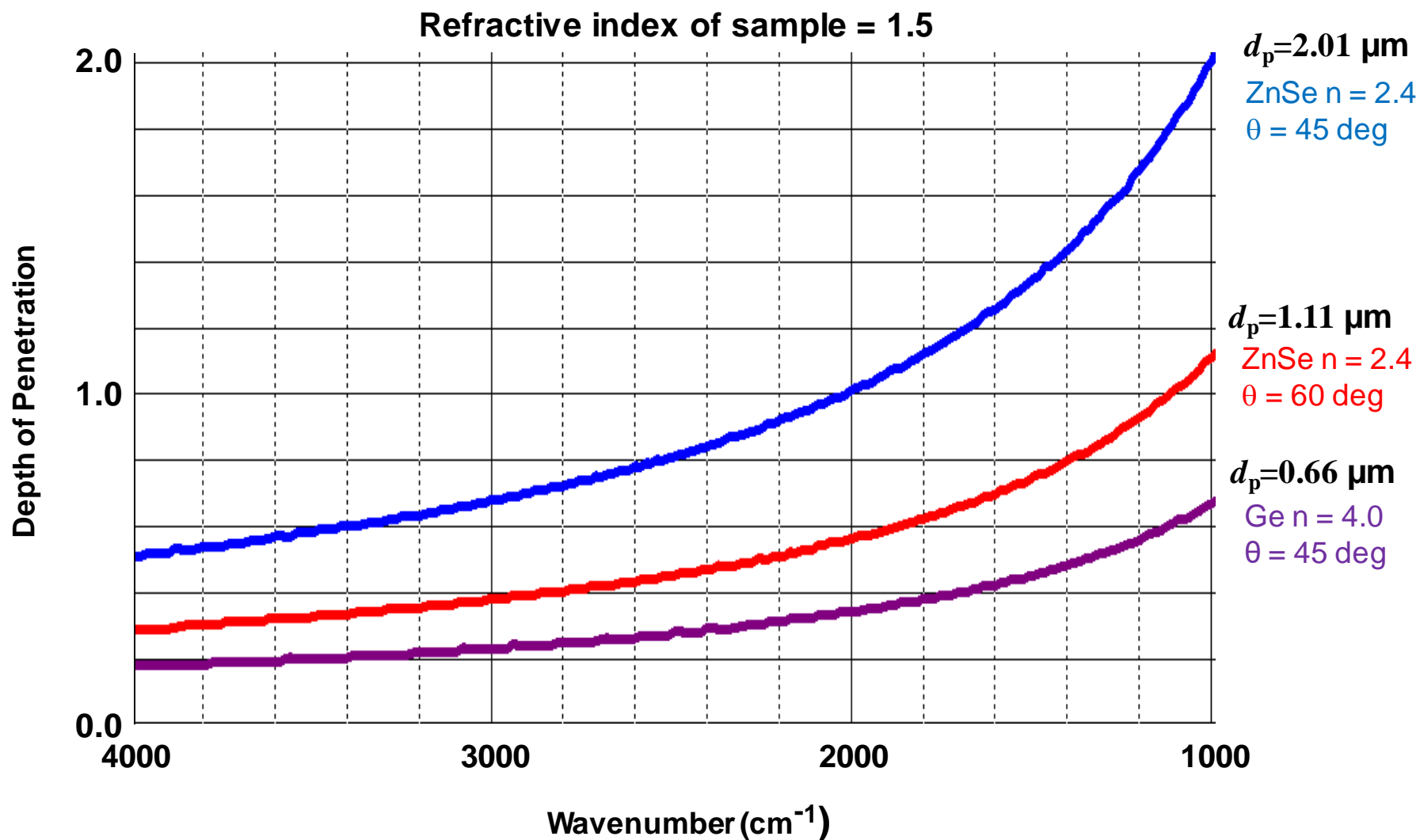

Failure Analysis via ATR


- Residue collected on filter membrane
- Analyzed by direct contact with Diamond ATR
 - Membrane + residue
 - Subtract membrane
 - Search

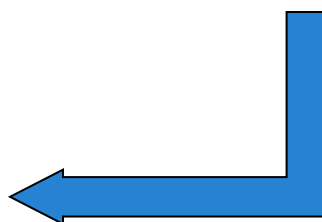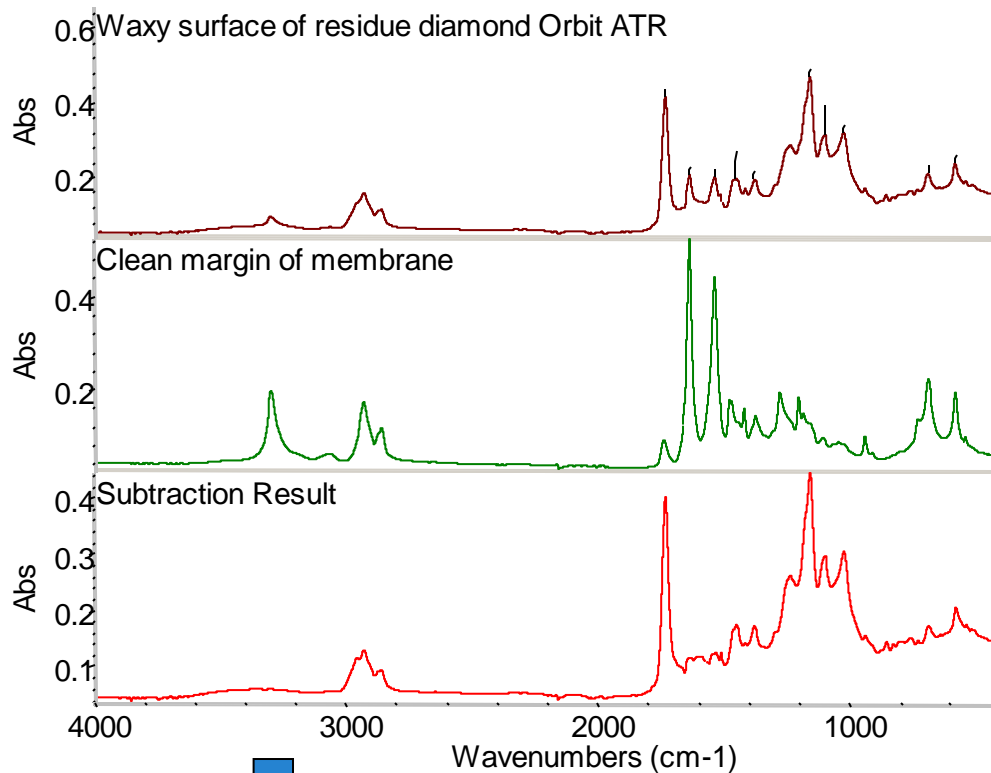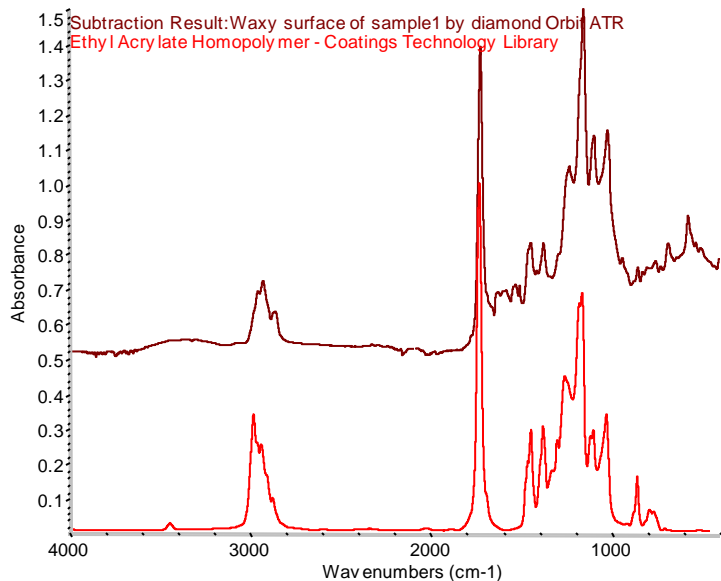

Ethyl Acrylate Homopolymer
Coatings Technology Library



Inorganic Residue on Turbine

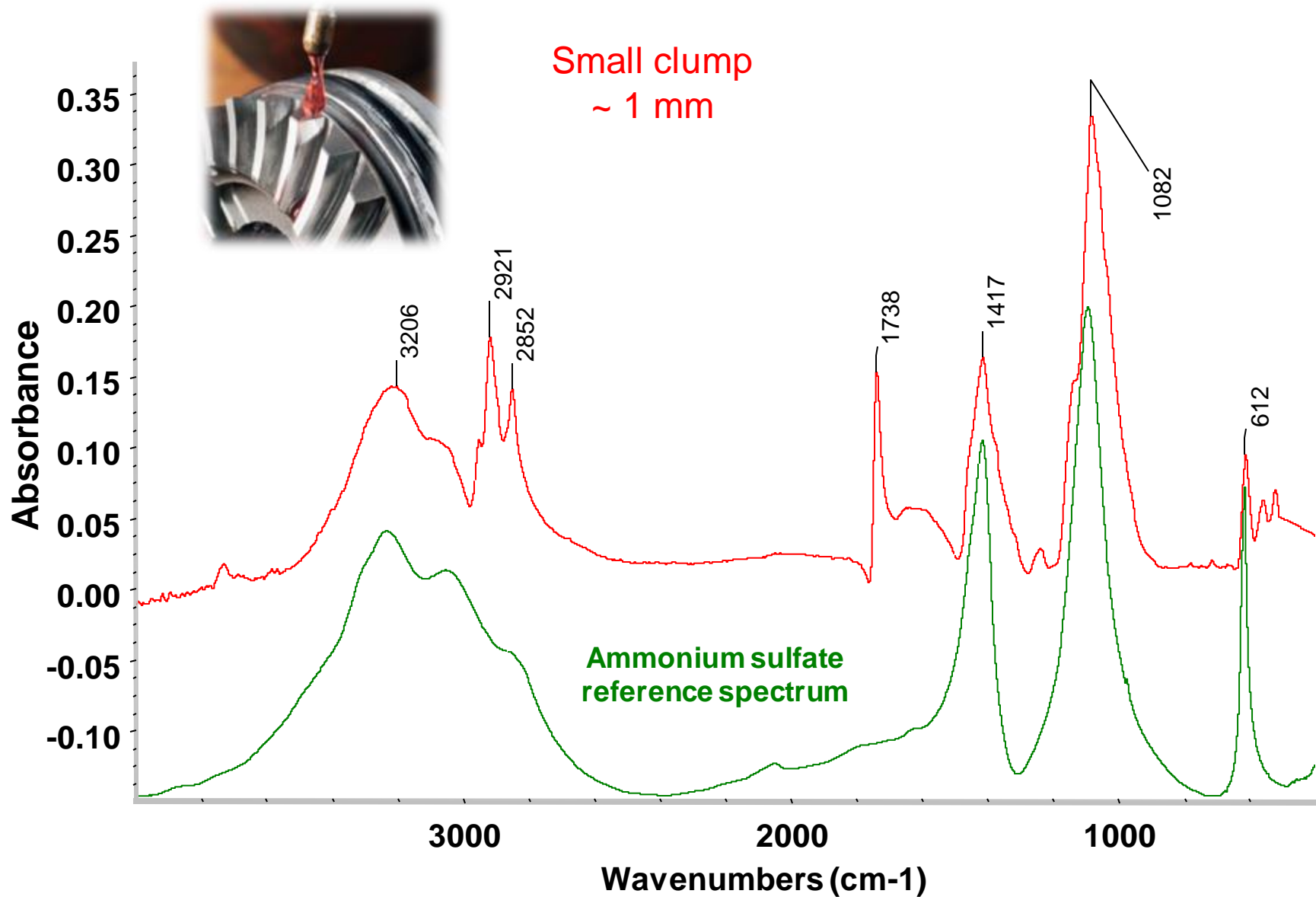


Heat Shrink Tubing

- Attempt to chemically identify and reverse engineer heat shrink tubing
 - Gardner Bender Polyolefin (HST-250)
- Since tubing appears to have a high loading of carbon black, analyze via germanium ATR as it has a lower depth of penetration



Initial Identification of Tubing


- Polyolefin tubing (red) library search result (blue) has strong match value
 - Polymer tubing C-H & C=O stretches appear to be highly correlated
- 3700 cm^{-1} stretch and low wavenumber region do not match well

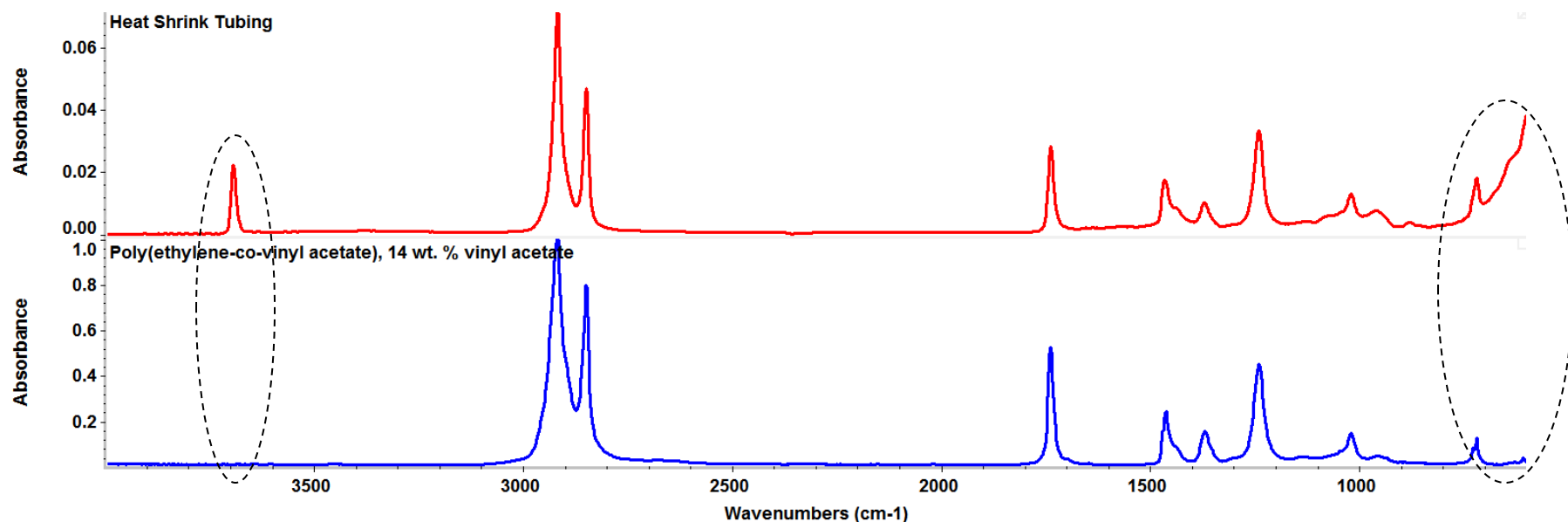

Index	Match	Compound Name	Library Name
1	17912	91.07 Poly(ethylene-co-vinyl acetate), 14 wt. % vinyl acetate	HR Aldrich FT-IR Collection Edition II
2	218	91.07 POLY(ETHYLENE-CO-VINYL ACETATE), 14 WT. % VINYL ACETATE	HR Aldrich Polymers
3	232	89.68 ELVAX 360	HR Industrial Coatings
4	119	89.58 ELVAX 420	HR Industrial Coatings


OMNIC Spectra Shows Multicomponent Match

- Peak at 3700 cm^{-1} and low wavenumber tail appears to be from an inorganic filler - magnesium hydroxide
- MgOH is used as a flame retardant

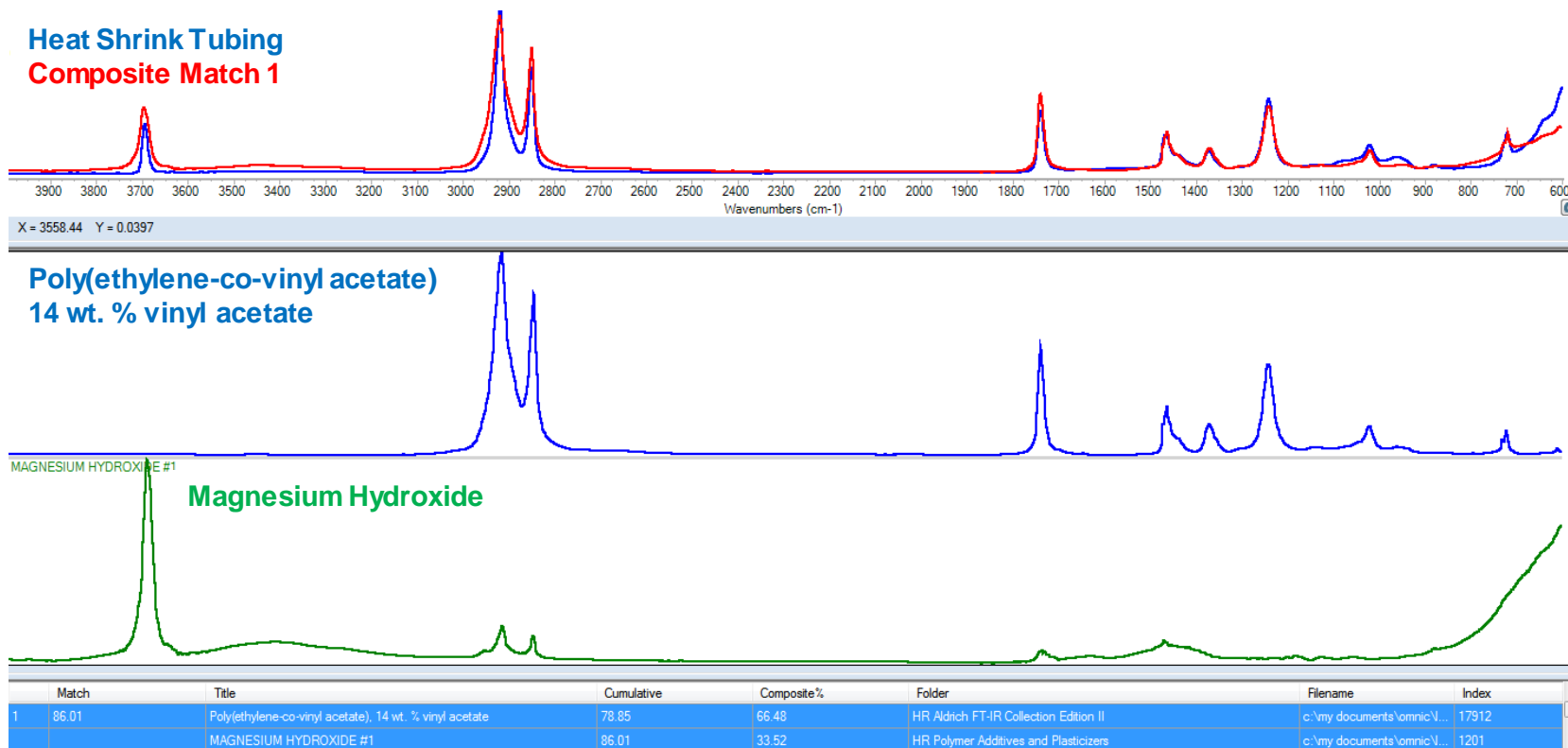


Outer Tubing

- Spectral changes from **inner tubing** is different from **outside** of the tubing
- Why would outside tubing have another chemical coating?
 - What is coating? And what is its purpose?

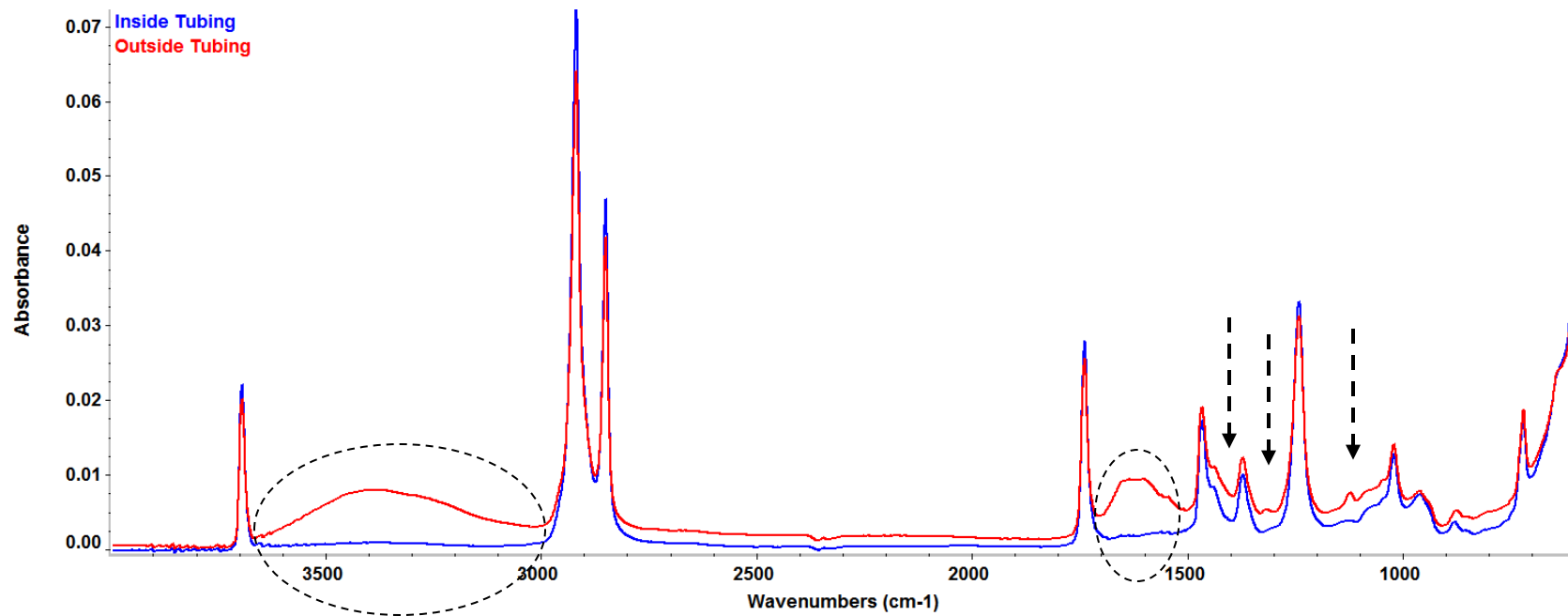



Identification of OD Coating

- Library search of outside coating shows match for Tyzor 131
 - Organic titanium complex made by DuPont as a cross-linking agent
- As polyolefin tube is heated – Tyzor 131 enhances crosslinking causing tube to shrink

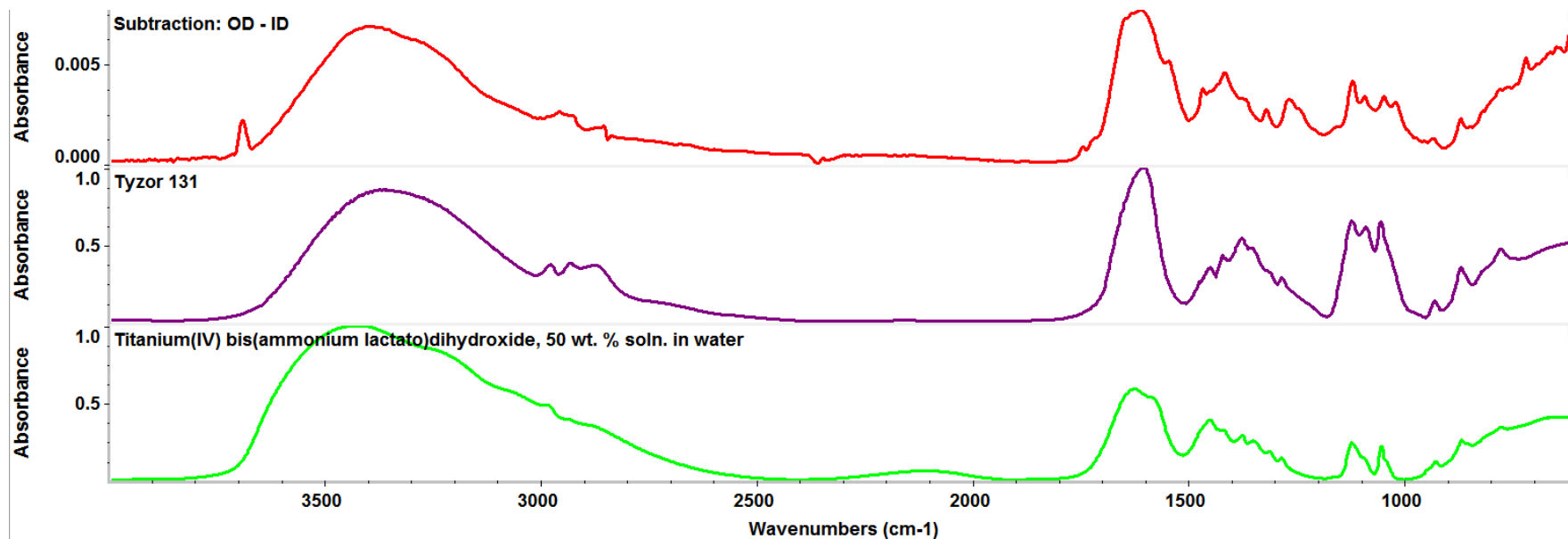

Index	Match	Compound Name	Library Name	
1	1671	59.90	Tyzor 131	HR Coatings Technology
2	17487	51.77	Titanium(IV) bis(ammonium lactato)dihydroxide, 50 wt. % soln	HR Aldrich FT-IR Collection Edition II
3	548	47.83	METHYL ALKYL IMIDAZOLINE SODIUM SALT	HR Polymer Additives and Plasticizers
4	573	47.28	Poly(terephthalovi oxamidrazone)+SrCO3	HR Hummel Polymer and Additives

TGA-IR: The Basics

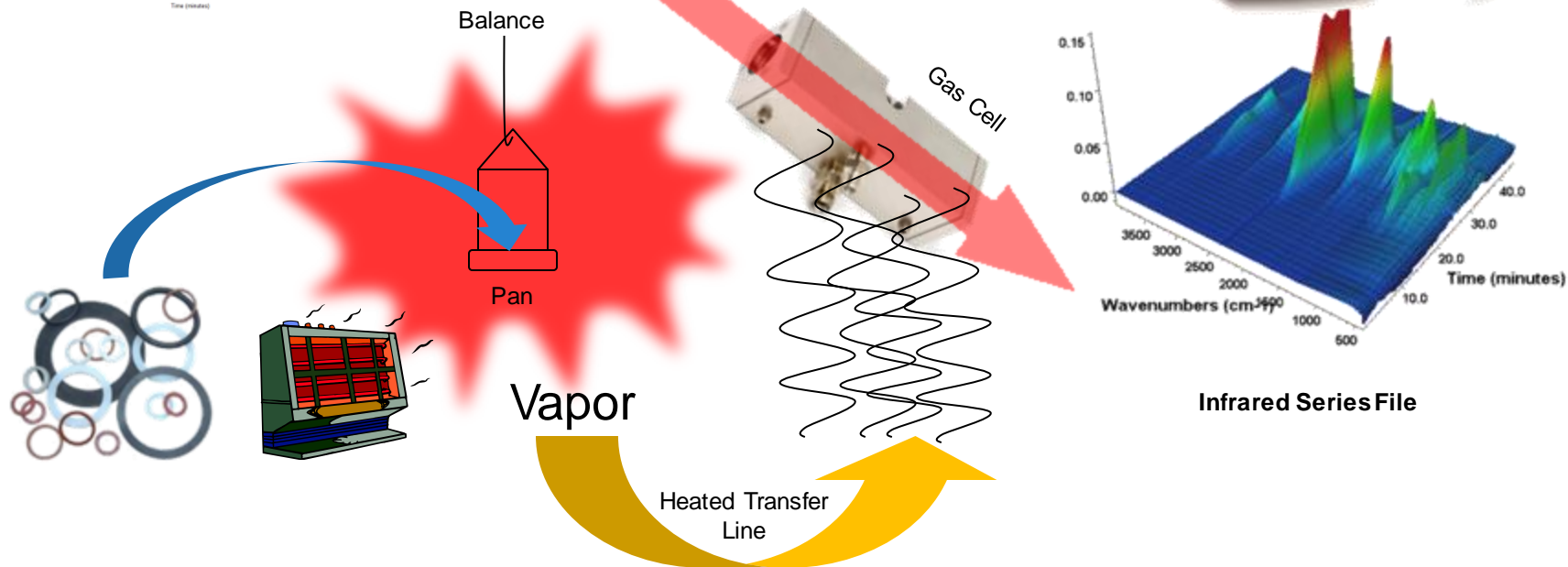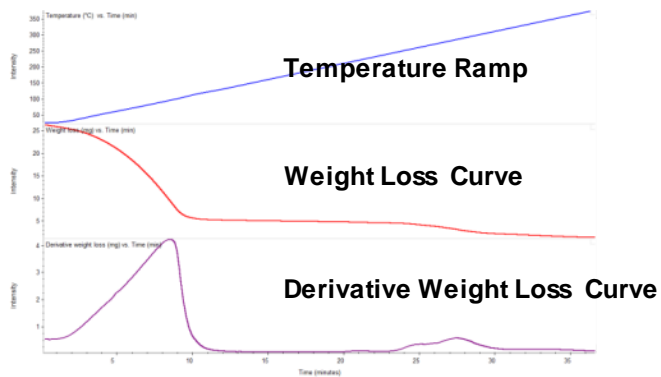

Quantitative: T "G" A



Qualitative: FT-IR




- Deformulation and Failure Analysis
 - Rubber gaskets were found to be failing in the field
 - TGA analysis of failing gaskets shows incorrect formulation
 - Subtraction result shows chemical differences

IR of vapor at 67 min

Gram-Schmidt Plot

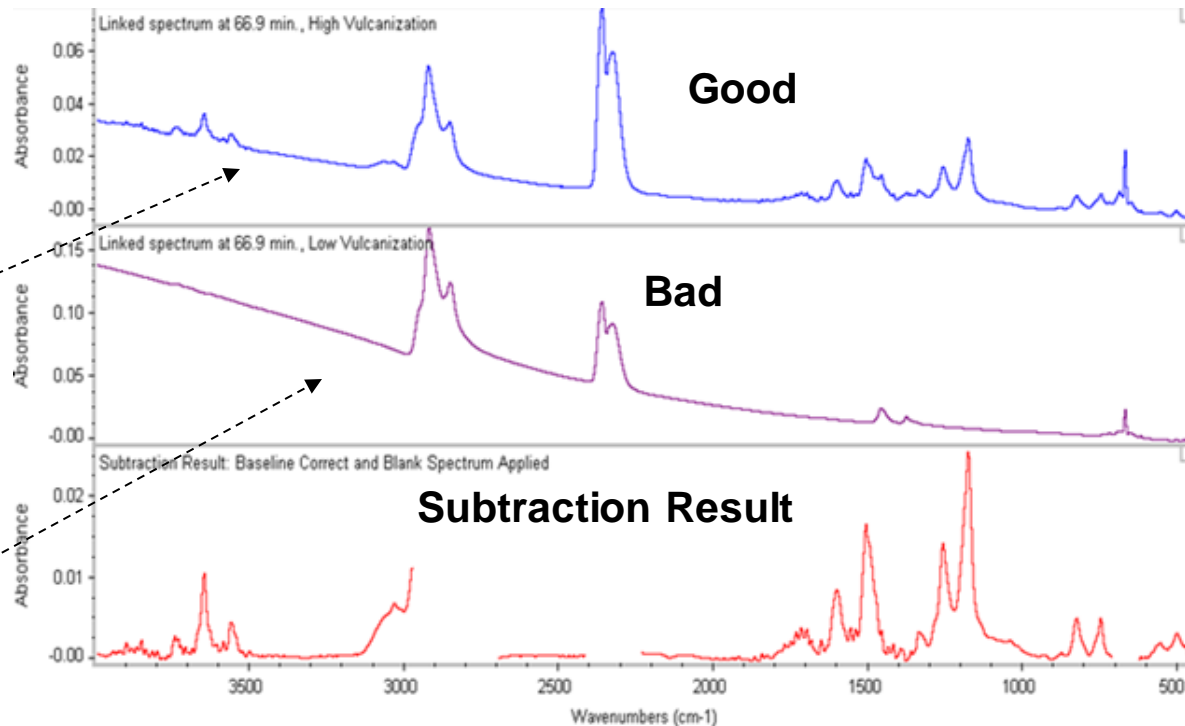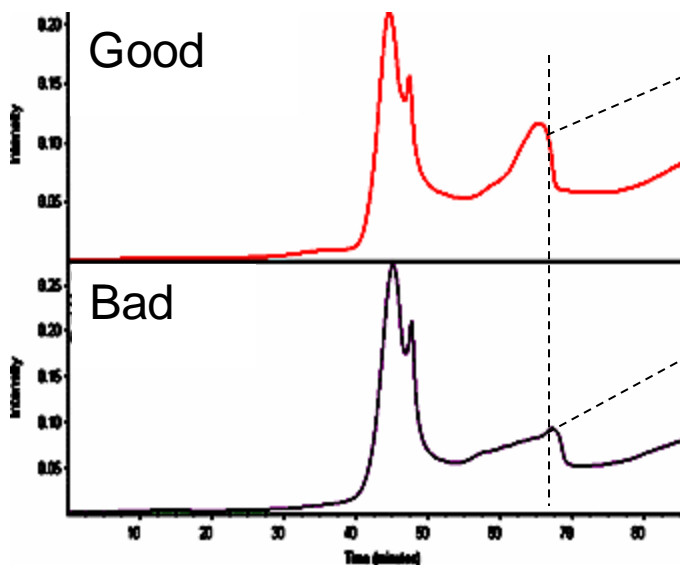


- Bad gasket does is missing bisphenol A
 - Chemical information to support early failure in field
- TGA coupled with IR allows for chemical evidence of failure

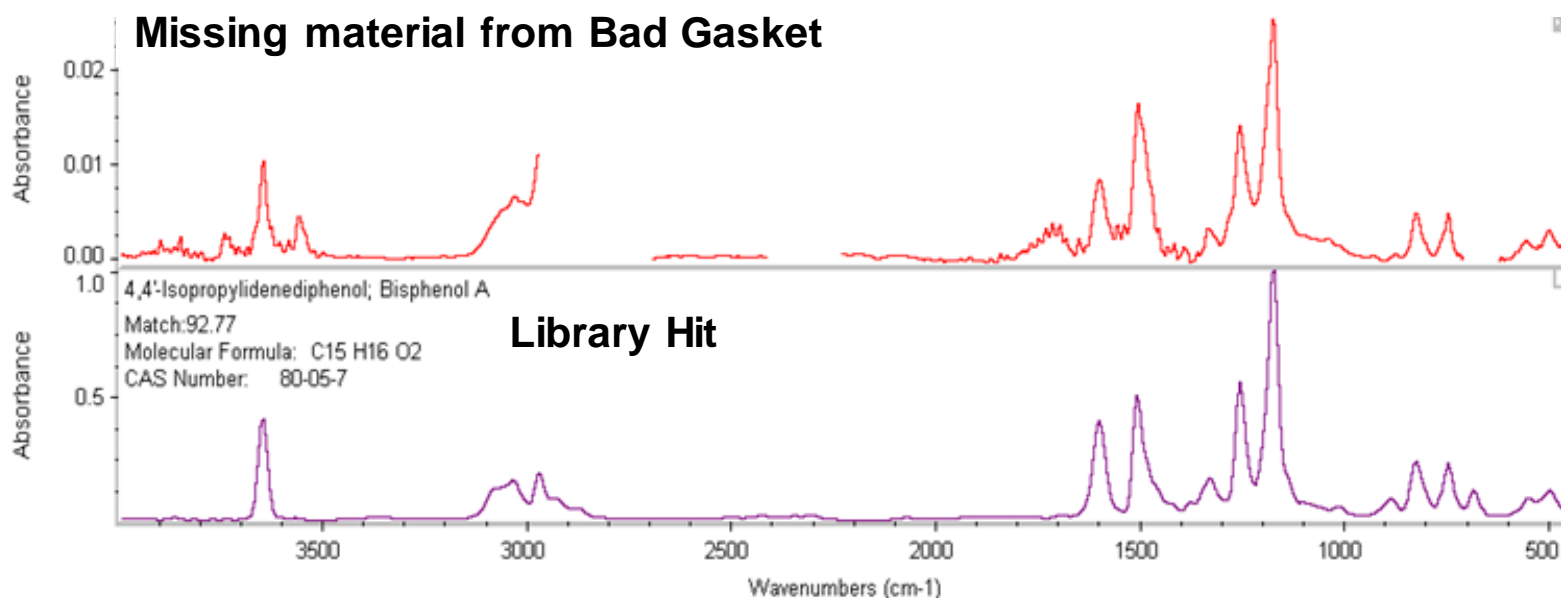


	Index	Match	Compound Name	Library Name	
	1	3269	92.77	4,4'-Isopropylidenediphenol; Bisphenol A	Nicolet/Aldrich Vapor Phase
	2	28	87.20	4,4'-Isopropylidenediphenol; Bisphenol A	Hazardous Chemicals - Vapor Phase Spectra
	3	5307	87.17	4,4'-Isopropylidenediphenol; Bisphenol A	Nicolet Vapor Phase
	4	407	87.17	4,4'-Isopropylidenediphenol; Bisphenol A	Nicolet TGA Vapor Phase
	5	5306	85.83	4-(2-Phenylisopropyl)phenol; 4-Cumylphenol	Nicolet Vapor Phase

Summary


- ATR was used to reverse engineer heat shrink tubing
- Both organic and inorganic components were identified

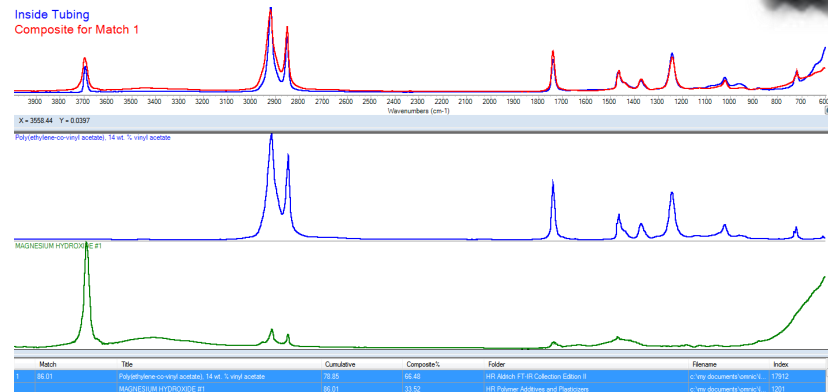

- TGA-IR aided in understanding chemical failure
- Chemical data coupled with physical data aids in understanding failure

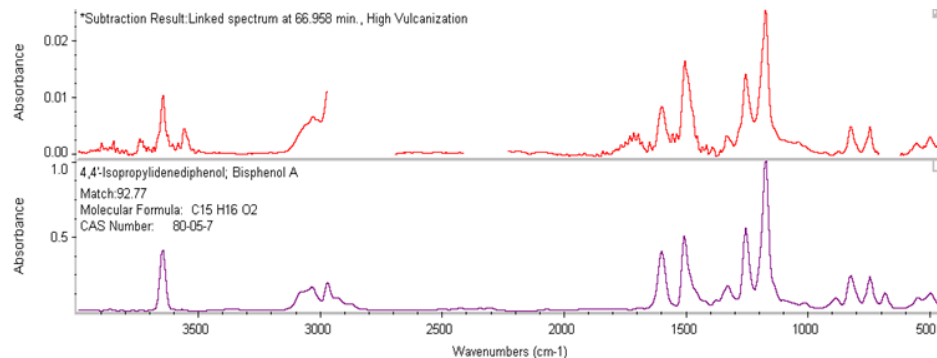

Thanks!

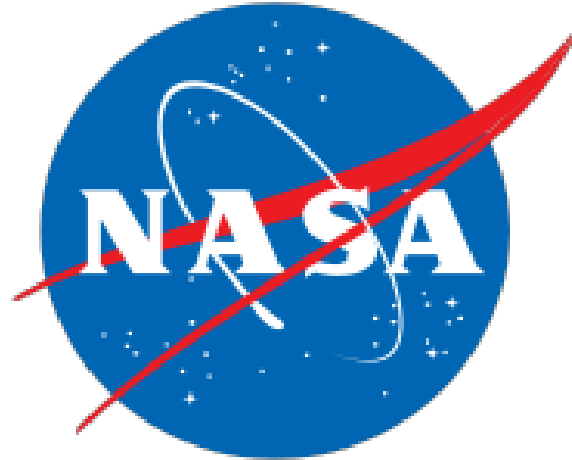

2017 NASA Contamination, Coatings, Materials, and Planetary Protection
Workshop

NASA Goddard Space Flight Center

Questions/Comments?